


AA / ICP / ICP-MS

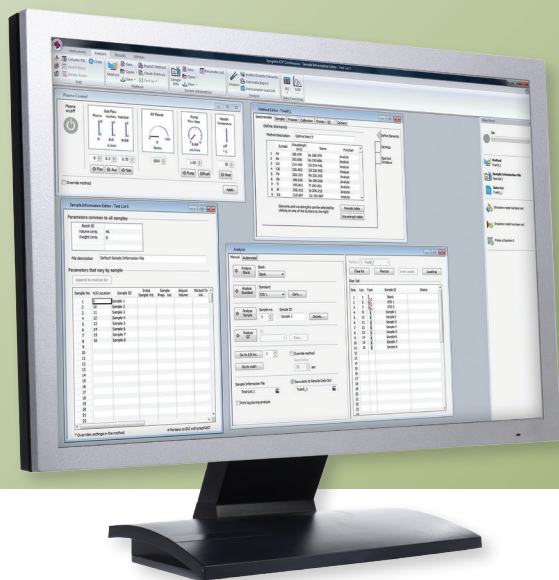

Key Capabilities

- Multiple user levels with login/password support through Microsoft® Windows®
- Method (and IEC or MSF files in the case of ICP) saved with the data in the results database
- Reprocessing doesn't change stored data
- Master event log records all significant actions performed by each user
- Files and data objects are automatically assigned version numbers and changes between versions are recorded

Syngistix Enhanced Security Software Speeds and Simplifies Regulatory Compliance

PerkinElmer's Syngistix™ Atomic Spectroscopy Software platform, the industry's first cross-platform instrument control solution, allows lab professionals to work seamlessly across multiple techniques, enhancing laboratory productivity. Plus, its unique icon-based design simplifies navigation and walks the user through every step of the analysis—from setting up to acquiring data to reporting results. Now an Enhanced Security™ version is available which meets the special needs of highly regulated labs such as those that must comply with the U.S. FDA's 21 CFR Part 11 regulations.

Standard Syngistix Package

Behind the intuitive interface of Syngistix lies a full suite of features designed to optimize data security and ensure compliance to help today's laboratories cope with the regulations mandated by government agencies or quality protocols.

- Syngistix leverages the powerful security features of Microsoft® Windows® to provide the protection your laboratory needs. Personnel can be divided into groups with each group assigned a level of permission. The "administrator" group can install and configure the software and perform any Syngistix task while other groups may be restricted from performing certain tasks. With Syngistix software, you define the groups and permissions assigned to each – passwords control access (Figure 1).

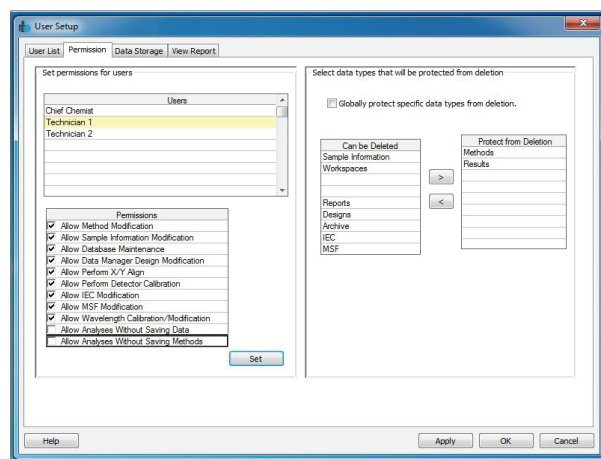


Figure 1. Syngistix user setup utility.

- Different users can access the same or different files such as method databases or sample information files as determined by the system administrator. Files can be located on the instrument controller or on other computers connected to the controller across the network.
- Syngistix “signs” its data. When data are saved from an analysis, a proprietary check sum is also computed and stored in the database – all transparent to the user. If anyone alters the data after it is saved, for example by altering the analysis time or changing the concentration value, this alteration is readily detected using a command in the Data Manager.
- Syngistix saves a copy of the method (and IEC or MSF files in the case of ICP) used to acquire the data along with the data in the results library. This provides an audit trail of the conditions used to perform the analyses.
- Reprocessing doesn’t change the data stored, but rather new data are written to the results library along with a notation stating that the data represent reprocessed data rather than original data.
- Syngistix software meets 21 CFR Part 11 requirements for a closed system.

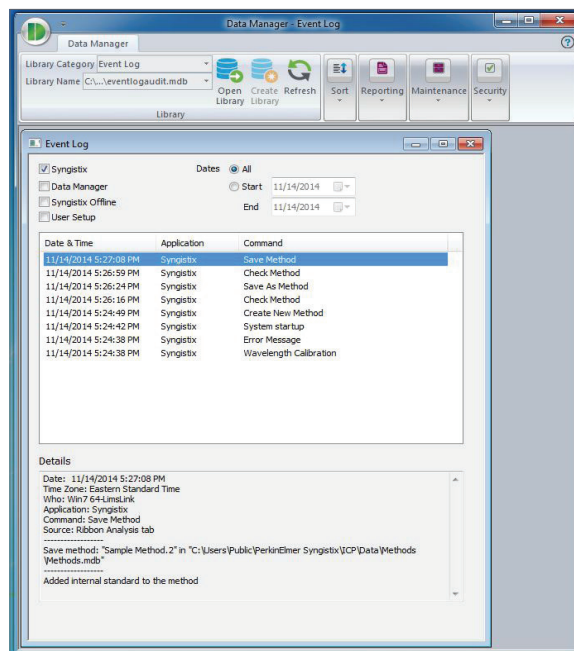


Figure 2. Viewer for the Master Event Log.

Enhanced Security Option

The Enhanced Security option adds the additional capabilities needed by highly regulated laboratories, such as those operating under the rules of 21 CFR Part 11. These additional capabilities supplement the features in the basic Syngistix package and include the following:

- The software maintains a Master Event Log that records all significant actions performed by the user. Each entry includes the date and time of the action, what was done, the name of the user, and, in many cases, the reason the action was performed. This event log can be viewed or printed using the Data Manager (Figure 2).
- The software automatically adds version numbers to all files and data sets, records the changes between versions, and automatically moves old versions to a history directory. Changes can be viewed using the Data Manager (Figure 3). The software restricts all users to choosing file names from a common pool. In this way, files are uniquely identified across the system, regardless of the analyst creating or using them.
- The software includes options to prevent analyses from being performed without saving data and to allow analyses only with saved methods. These restrictions ensure that a proper audit trail is maintained for all activities.

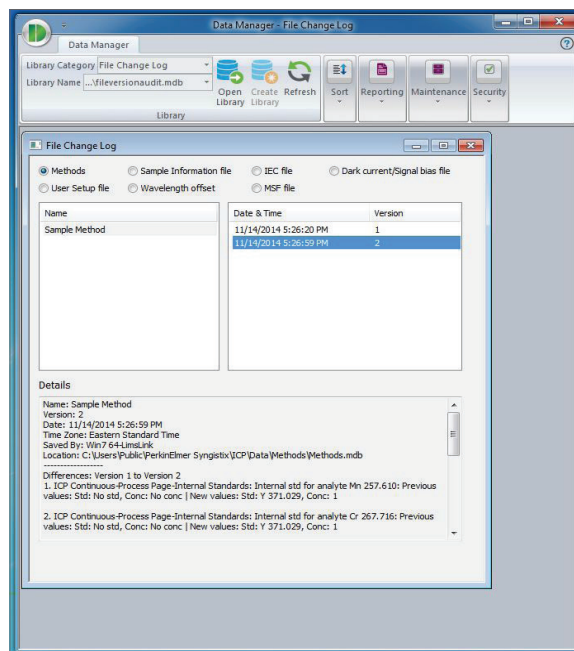


Figure 3. Viewer for the File Change Log where differences between versions of files and data objects are displayed.